

USER GUIDANCE

SUITE[®] TOUCHLESS FRAME WITH REPLACEABLE NYLON HOOKING STRIPS

GRADE C AND D NON-STERILE MANUFACTURING:

- 1- Open bag with clean and individually packaged Suite[®] Frame with attached nylon strips
- 2- Clean and/or disinfect environment with Suite[®] Touchless Mops
- 3- Rinse Suite[®] Frame fully after each use
- 4- Repeat steps above for each use
- 5- Replace strips in frame as needed, usually 3 months depending on use
- 6- Replace Suite[®] Frame as needed, usually 1 to 2 years depending on usage



GRADE A AND B STERILE MANUFACTURING:

- 1- Autoclave Suite[®] Frame prior to use
- 2- Place pre-sterilized nylon hooking strips into frame, put end cap on frame
- 3- Clean and/or disinfect sterile environment with Sterile Suite[®] Touchless Mops
- 4- Rinse Suite[®] Frame fully after each use, discard strips
- 5- Repeat steps above for each use
- 6- Replace Suite[®] Frame as needed, usually 6 months to 2 years depending on usage

