



DPTE-BetaBag®
Standard Range PU Bag
Gamma and X-Ray Sterilization Range

Product Sheet – FP 213 | 006

Product overview

Application

The DPTE-BetaBag® is a single use solution for safe, fast, contamination free transfer in and out of an isolator or filling line.

The DPTE-BetaBag® with a PU film bag is mainly used for waste disposal, but in some cases inlets compatible with gamma and X-Ray sterilization, are possible.



Key features

DPTE-BetaBag® Part Number	DPTE® BETA Type	Useful Volume	Bag size mm (LxW)	Drawing number	Sterile	
					YES	NO
6007001617	105	30 L	930 x 400	4090-13-01	X	
6007001925	190	10 L	550 x 310	4090-23-28	X	
6007001926	190	10 L	550 x 310	4090-23-28		X
6007001989	190	30 L	1000 x 400	4090-23-26	X	
6007001990	190	30 L	1000 x 400	4090-23-26		X
6007002098	190	100 L	1150 x 740	4090-23-27	X	
6007002097	190	100 L	1150 x 740	4090-23-27		X
6007001829	190	150 L	1400 x 740	4090-23-24	X	
6007001624	190	150 L	1400 x 740	4090-23-24		X
6007002345	270	35 L	730 x 558	4090-33-10	X	
6007002344	270	35 L	730 x 558	4090-33-10		X
6007001105	270	100 L	1150 x 815	4090-33-09	X	
6007002346	270	100 L	1150 x 815	4090-33-09		X
6007002348	270	150 L	1300 x 740	4090-33-08	X	
6007002347	270	150 L	1300 x 740	4090-33-08		X

Product description

Main components and materials

DPTE-BetaBag® = DPTE® Beta part + Bag

- **DPTE® Beta part**
 - DPTE® Beta flange: HDPE material
 - DPTE® Beta door: HDPE material
 - J3L Silicone gasket.
 - Protective cover (J3L gasket protection against damage and dirt) made with:
 - PVC material for Ø 105 and Ø 270.
 - PC material for Ø 190.
- **Bag**
 - Body: Polyurethane film 200 µm.
 - No inner sleeve.
 - Cylindrical tight sealing onto DPTE® Beta part.

Product specificities

- Sterilization indicator located onto the first packing and on the cardboard box for sterile products.
- Single use DPTE-BetaBag®.
- Multi-connection on DPTE® system (qualified to be connected/disconnected 5 times).
- For non sterile product:
 - Compatible for gamma and X-Ray sterilization.
 - Designed to be gamma or X-Ray sterilized once.

Loading

- Loading of the DPTE-BetaBag® has to be performed through the DPTE® Beta Part connected onto a DPTE® Alpha Part.
- Maximum recommended load: 10 kg (depending on type of components).
- We suggest using a trolley when handling loaded bags.

Sterility

Sterile products, packed according the table “packaging” have been qualified according to standards ISO 11137-1 and 2 and ISO 11737-2 for a gamma sterilization cycle between 25 and 50 kGrays.

Properties

DPTE-BetaBag® PROPERTIES

PROPERTIES	TEST METHOD	UNITS	ACCEPTANCE CRITERIA
Leaktightness <ul style="list-style-type: none">• DPTE® Beta Part 105• DPTE® Beta Part 190• DPTE® Beta Part 270• DPTE-BetaBag®	<p>Pressure differential at P= -4000 Pa</p> <p>Pressure differential at P=1000 Pa</p>	<p>Pa/min</p> <p>Pa/min</p>	<p>According to Getinge specification</p>
Seal strength	ASTM F88	N/15 mm	≥ 15

PU FILM PROPERTIES

Refer to FT 096 – PU film data sheet

Manufacturing process

Cleanroom classification is according to ISO 14644-1 standard.

- Shaping of the liner in ISO 7 cleanroom.
- Tight assembly of the bag onto the DPTE® Beta flange in ISO 7 cleanroom.
- Assembly of the J3L gasket and DPTE® Beta door in ISO 7 cleanroom.
- Packaging in double bag in ISO 7 cleanroom.

Routine tests

- **Tests performed 100 % in ISO 7 cleanroom.**
 - Pressure differential test of the DPTE® Beta Part.
 - Pressure differential test of the DPTE-BetaBag®
- **Sampling test** (based on Standard NF ISO 2859-1) performed in ISO 7 cleanroom.
 - Dimensional and visual check.

Traceability

- Each DPTE-BetaBag® is identified by a batch number.
- Each batch number is associated with a manufacturing batch record.

Safety features

- The design of the DPTE-BetaBag® ensures that it cannot be connected to the DPTE® Alpha part if the DPTE® Beta door is not properly closed.

Product compliance

- **DPTE® Beta flange, door**
 - Meets the FDA 21 CFR 177.1520 regulation
 - BSE/TSE: EMEA 410/01 compliant
- **PU film**
 - Meets the FDA 21 CFR 177.2600
- **J3L silicone gasket Ø 105**
 - Meets the FDA 21 CFR 177.2600 regulation
- **J3L silicone gasket Ø 190**
 - Meets the biocompatibility USP <88> Class VI
 - Meets the FDA 21 CFR 177.2600 regulation
 - BSE/TSE: the material contains no substances of animal origin.
- **J3L silicone gasket Ø 270**
 - Meets the biocompatibility USP Class VI
 - Meets the FDA 21 CFR 177.2600 regulation

Documentation

A statement of conformity is delivered with each manufacturing batch.

A gamma irradiation treatment certificate is delivered with each sterile manufacturing batch.

Use

Service conditions

- Operating temperature for loading and unloading: between 16 and 24 °C.
- Gamma and X-Ray sterilization dose validated between 25 and 50 kGrays.

Limits of resistance

Pressure: Maximum + 1000 Pa (relative pressure).

Guarantee and shelf life

Under normal conditions of use and storage.

- DPTE-BetaBag® Ø 190: 30-month shelf life after manufacturing date.
- DPTE-BetaBag® Ø 105 and Ø 270: 24-month shelf life after manufacturing date.

Shelf life validation ongoing for X-Ray range.

Accessories for use (handles)

- DPTE-BetaBag® 105: manual handling (no tool required).
- DPTE-BetaBag® 190 and 270: Set of 2 handles.

P/N 35502 C



Compatible type of DPTE® Alpha

DPTE®-XS with safety interlocks (FP 211).

Packaging & Storage

Packaging

DPTE BetaBag® type	105	190	270
Double over packaging	By 4 units	By 3 units	By 1 unit
Cardboard box	6 packs of 4 units	3 packs of 3 units	6 packs of 1 unit
Cardboard dimensions (mm)	780x420x360	780x420x360	550x385x570
Bag per box	24 units	9 units	6 units
Boxes per pallet	15 boxes		
Pallet size over all (mm)	1200x1000x1950		
Bag per pallet	360 units	135 units	90 units

Packaging identification

- Item designation
- Item reference number
- Batch number
- Quantity

Storage conditions

- Store the DPTE-BetaBag® in its original packaging between 16 and 24 °C and at relative humidity < 70 %.
- Protect the DPTE-BetaBag® against the effects of strong light (sun, electric lighting) and strong heat (radiators, electrical heaters).
- Use the DPTE-BetaBag® on first-in, first-out basis.

Glossary

DPTE®

“Double Porte de Transfert Etanche” Double door for leaktight transfer

DPTE®-XS

eXtra Secure Double door for leaktight transfer

HDPE

High Density PolyEthylene

J3L

Triple effect lip seal

PC

PolyCarbonate

PVC

PolyVinyl Chloride



With a firm belief that every person and community should have access to the best possible care, Getinge provides hospitals and life science institutions with products and solutions aiming to improve clinical results and optimize workflows. The offering includes products and solutions for intensive care, cardiovascular procedures, operating rooms, sterile reprocessing and life science. Getinge employs over 10,000 people worldwide and the products are sold in more than 135 countries.

Manufacturer:

Getinge Life Science France – Vendôme site

1 rue du Comté de Donegal – 41102 Vendôme cedex, France
Phone: +33 [0] 254 734 747 / Fax: +33 [0] 254 734 748

www.getinge.com